

Azura 2x HS Taq Mix

Catalog No.	Pack Size and Concentration	Components and Volume
AZ-1602	200 x 50 µl reactions, 2x	2x HS Taq Mix - 4 x 1.25 mL
AZ-1610	1000 x 50 µl reactions, 2x	2x HS Taq Mix - 20 x 1.25 mL
AZ-1620	2000 x 50 µl reactions, 2x	2x HS Taq Mix - 10 x 5 mL

Description

Azura 2x HS Taq Mix is a ready-to-use single tube hot-start PCR mix optimized for convenience and reduced pipetting steps. The highly advanced Azura 2x HS Taq Mix chemistry allows fast cycling conditions to be used with greater efficiency, reproducibility and throughput.

- New-generation PCR formulation including enhancers for maximum PCR efficiency, sensitivity and reaction speed.
- Robust Hot-Start amplification across a wide range of templates and sequence complexity including GC-rich DNA.
- Supplied as a ready-to-use 2x master-mix for maximum convenience and minimal liquid handling.
- High-yields with amplicons up to 5 kb with standard or fast cycling.

Storage

Azura 2x HS Taq Mix is shipped on blue ice and should be stored at –20°C upon receipt. Excessive freeze/thawing should be avoided.

Important Guidelines

Azura 2x HS Taq Mix: The 2x mix is comprised of Azura HS Taq DNA Polymerase, 2 mM dNTPs, 6 mM MgCl₂, and PCR enhancers for maximum efficiency, sensitivity and success with difficult amplicons. We do not suggest the use of additional PCR enhancers.

Template: For complex genomic DNA, we suggest the use of 5 ng – 500 ng per reaction; For cDNA or plasmid DNA, please use < 100 ng per reaction.

Primers: Primers should have a predicted melting temperature of around 60°C, using default Primer 3 settings (<http://frodo.wi.mit.edu/primer3/>). The final primer concentration in the reaction should be between 0.2 µM and 0.6 µM.

Annealing: We strongly recommend performing a temperature gradient to determine the optimal annealing temperature. Alternatively, 58°C can be used as a starting point. The optimal annealing temperature is usually 2-5 °C below the lower T_m of the pair. Depending on the reaction, the annealing time can also be reduced to 5 seconds.

Extension: Optimal extension is achieved at 72°C. The optimal extension time is dependent on amplicon length and complexity. For low complexity template such as plasmid DNA, an extension time of 15 seconds is sufficient for amplicons under 1 kb. For amplification from eukaryotic genomic DNA or cDNA, 30 seconds per kb is recommended.

Reaction setup: Allow Azura 2x HS Taq Mix to thaw and mix well by inversion. Centrifuge prior to use.

1. Prepare a PCR master mix based on following table:

Component	50 µl Reaction	Final Concentration/Notes
Azura 2x HS Taq Mix	25 µl	1x
Forward Primer (10 µM)	2.0 µl	400 nM
Reverse Primer (10 µM)	2.0 µl	400 nM
Template DNA	<100 ng cDNA, <500 ng genomic DNA	variable
PCR-grade water	Up to 50 µl final volume	

* For alternative total reaction volumes (eg. 20 µl), scale all components proportionally to maintain final concentrations.

2. PCR cycling:

Cycles	Temperature & Time	Notes
1	95°C, 2 minutes	Initial Denaturation, enzyme activation
25 - 40	95°C, 15 seconds 60°C, 15 seconds 72°C, 30 seconds per kb	Denaturation Annealing* (determined by user) Extension*

* See Important Guidelines.

Quality Control

Azura 2x HS Taq Mix is tested extensively for robust activity, processivity, efficiency, heat activation, sensitivity, absence of nuclease contamination and absence of nucleic acid contamination. Azura 2x HS Taq Mix is manufactured under a comprehensive quality management system, following ISO 9001:2015 standards.

Limitations of Use

This product is intended for research purposes only and is not intended for any animal or human therapeutic use.

Technical Support

For Trouble-shooting and Technical Guidance, please contact us at tech@azuragenomics.com and provide PCR reaction conditions, cycling parameters, amplicon size, and screen grabs (gel images) if possible.

Azura Genomics guarantees the performance of all products in the manner described in our product literature. The purchaser must determine the suitability of the product for its particular use. Azura Genomics shall not be liable for any direct, indirect, consequential or incidental damages arising from the use, the result of use or the liability to use this product.